




# 26400 WATER SATURATED TOLUENE VIAL DIRECTIONS FOR USE

## Step One

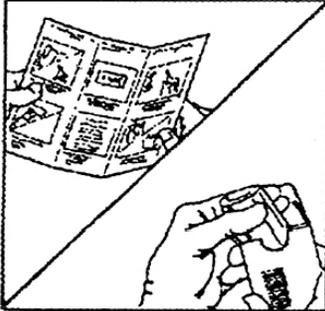


- Read and follow all directions carefully and confirm procedures with your supervisor.
- Wear protective eyewear, hand protection and safety clothing.
- Use in a well ventilated area.
- Break carton label seal and remove vial from the sealed bubble bag.

## Step Four




- Cut the shrink seal and pour the saturated Toluene as recommended in API MPMS Chapter 10.3, 10.4 and 18.1 into a volumetrically verified centrifuge tube.

## Step Two

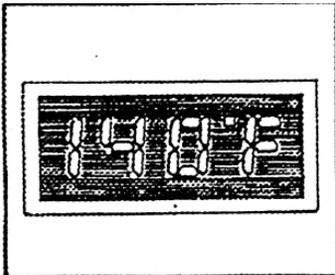

- Heat Vial to 140F (60C) in Oven or Heated centrifuge. (160F if using that method).
- Remove heated vial after 30 minutes.

## Step Five

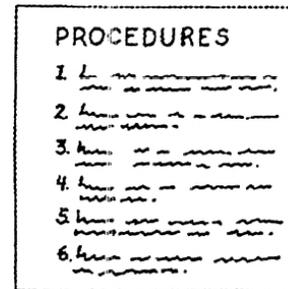


- Follow your test methods or your supervisor instructions.

## Step Three



- Shake Toluene vial until all free water is in suspension

## Step Six

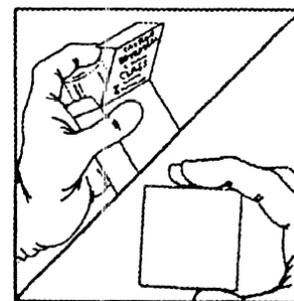

- Place empty vial and cap into the packaging and dispose of properly